


Upload process lists to spectrometer

Process Lists cannot be placed manually on the spectrometer. Even though they are in the expected location they will not appear in the “Process Lists File Browser” in Delta as the displayed listing is from a database, not from the physical location of files.


Disconnect from the Spectrometer before the following operation is performed.

Navigate to the location of the process lists – “proton_autophase_subset.list”,
“carbon_autophase_subset.list”.


Select “proton_autophase_subset.list”

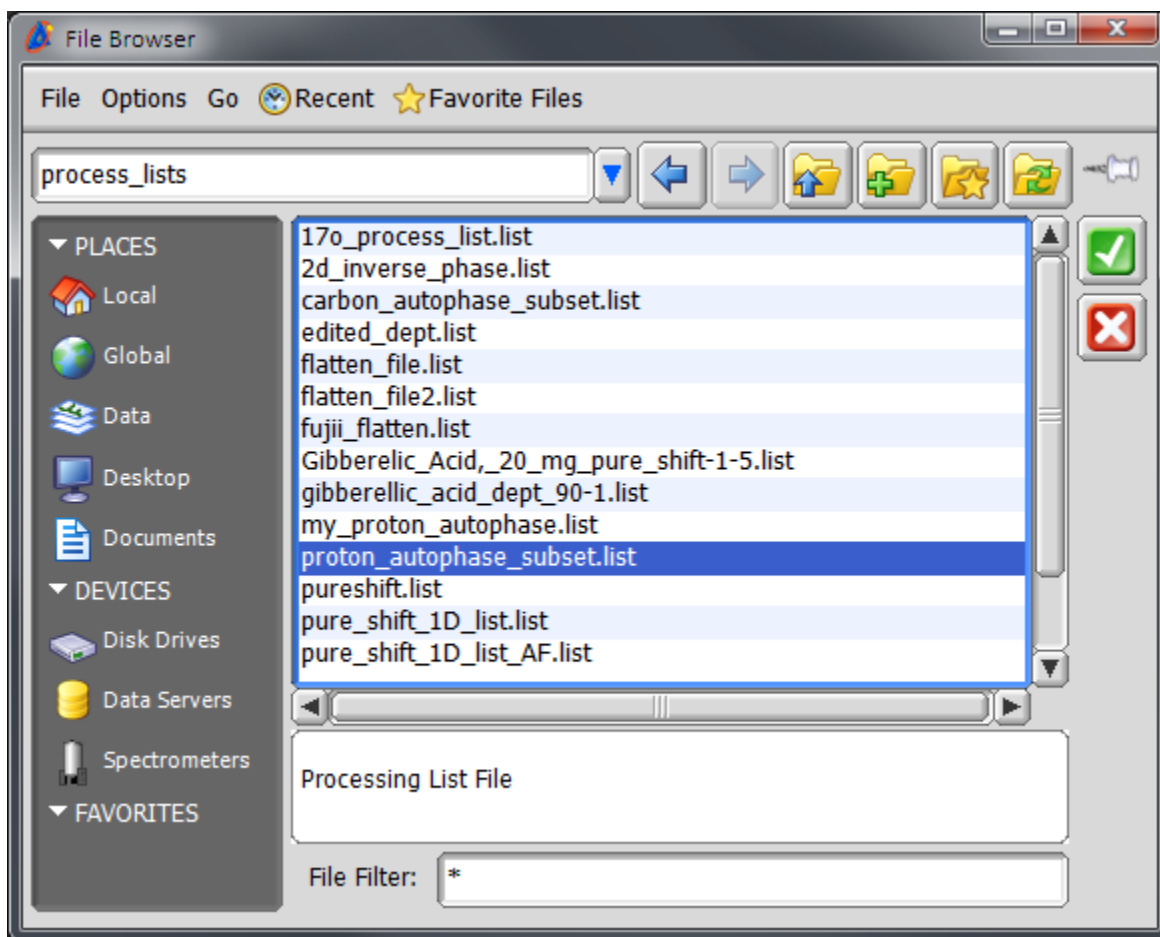


From the “File” menu select “Upload File”

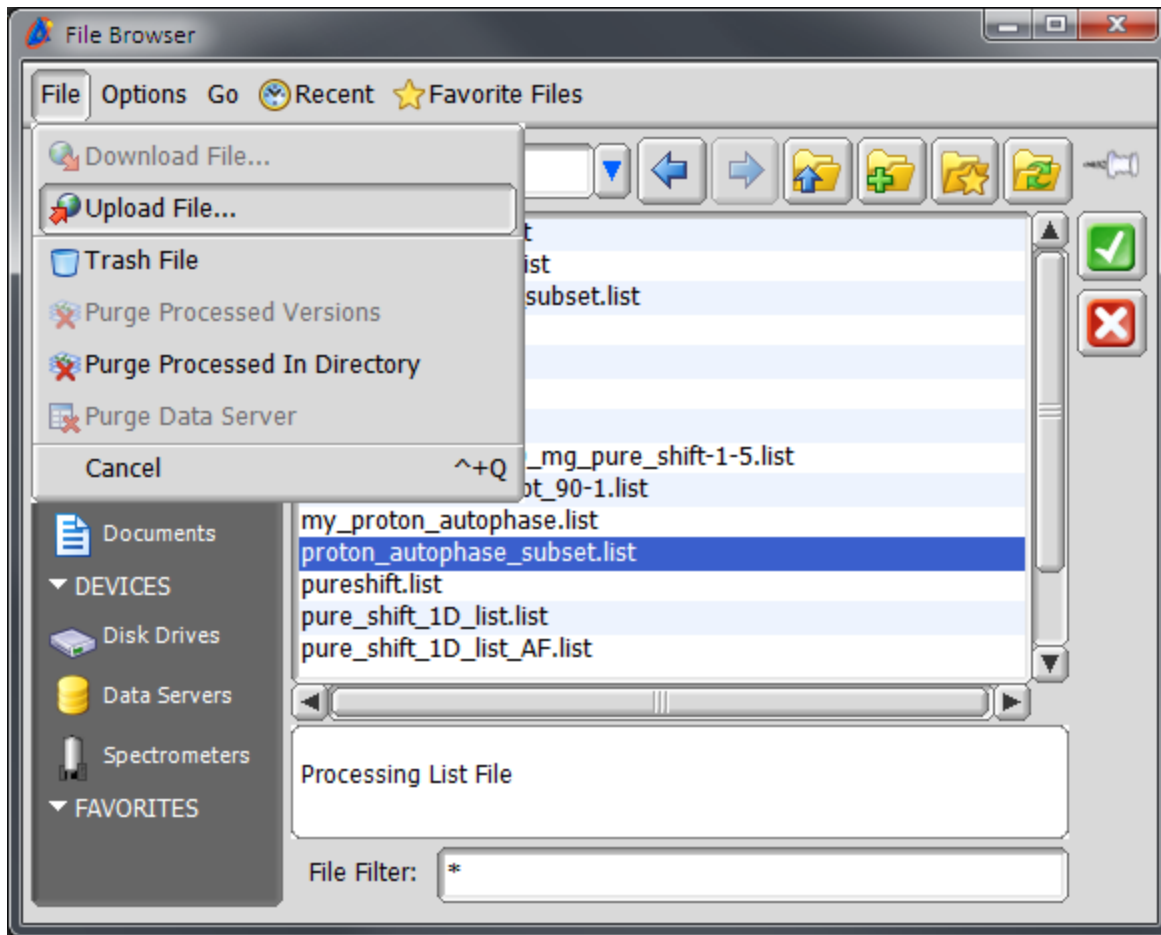

A new window displaying available spectrometers will appear. Select the spectrometer to which the file needs to be uploaded to.

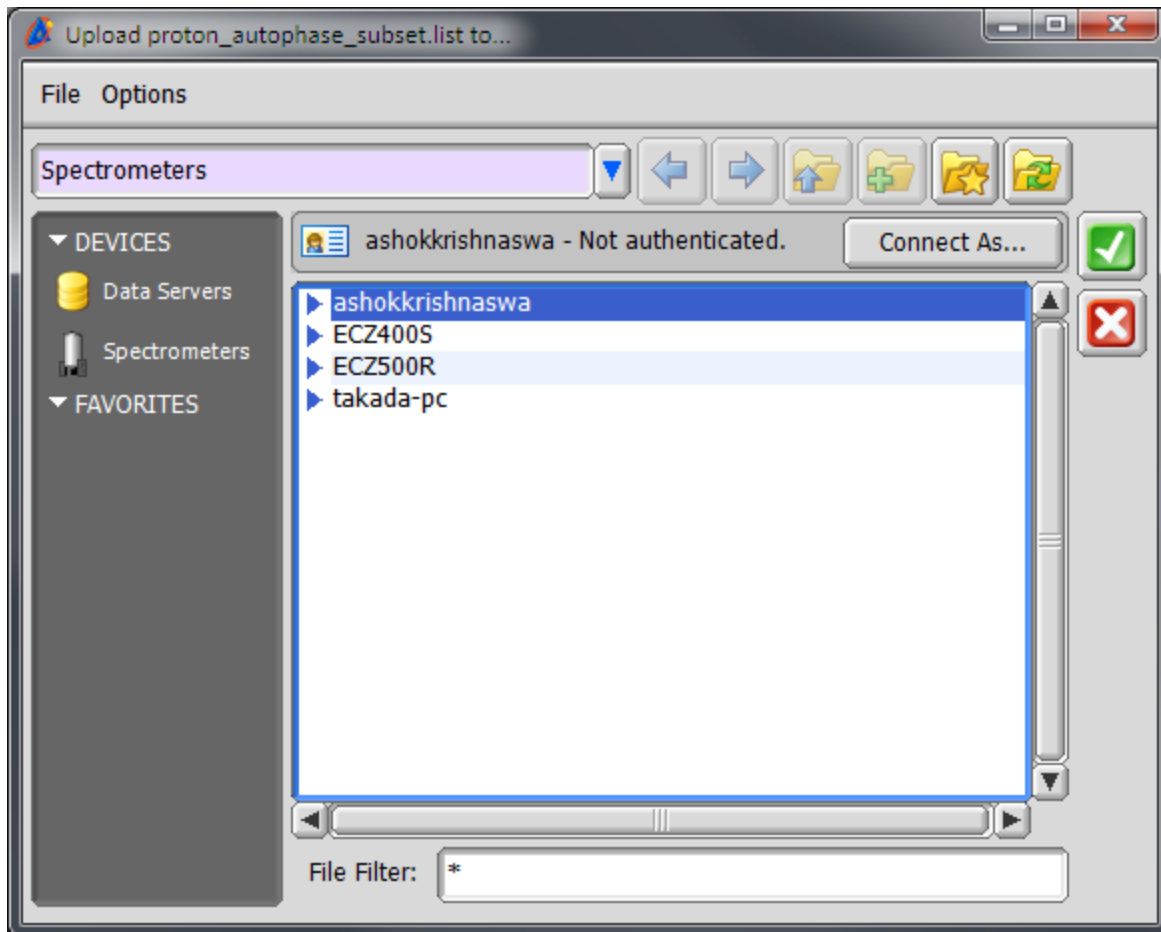

Click on "Connect As..."

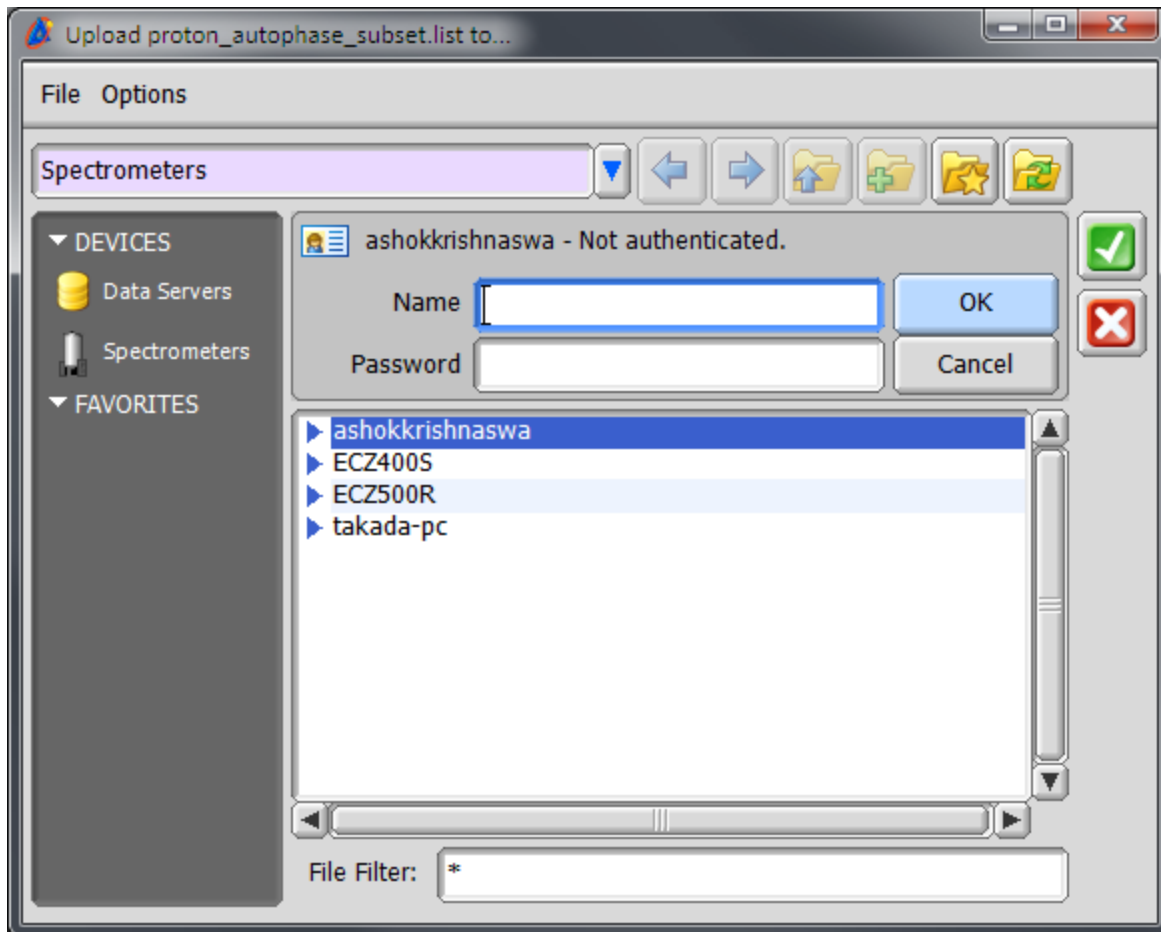


Enter


Name: console

Password: console


Hit the "OK" button.


After authentication double click on the spectrometer to show sub folders.

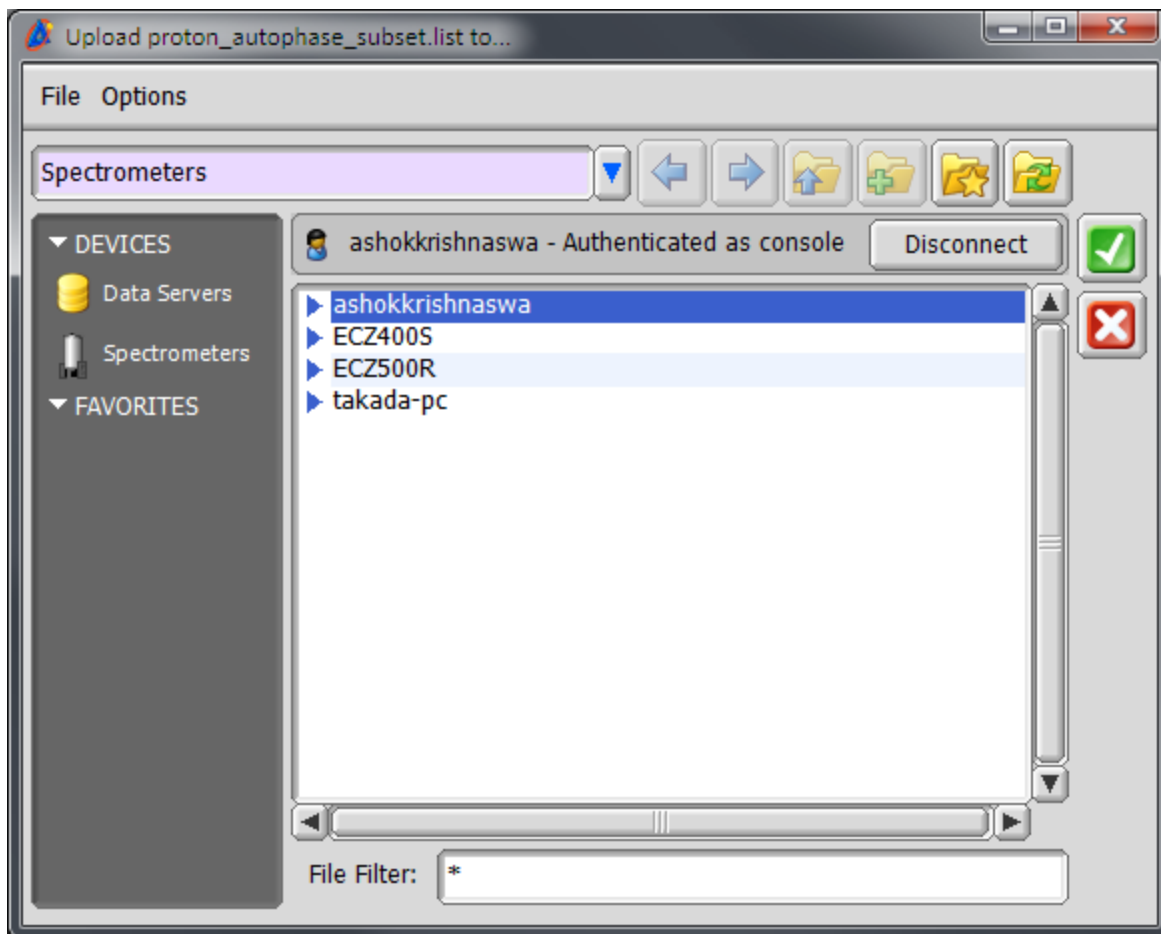

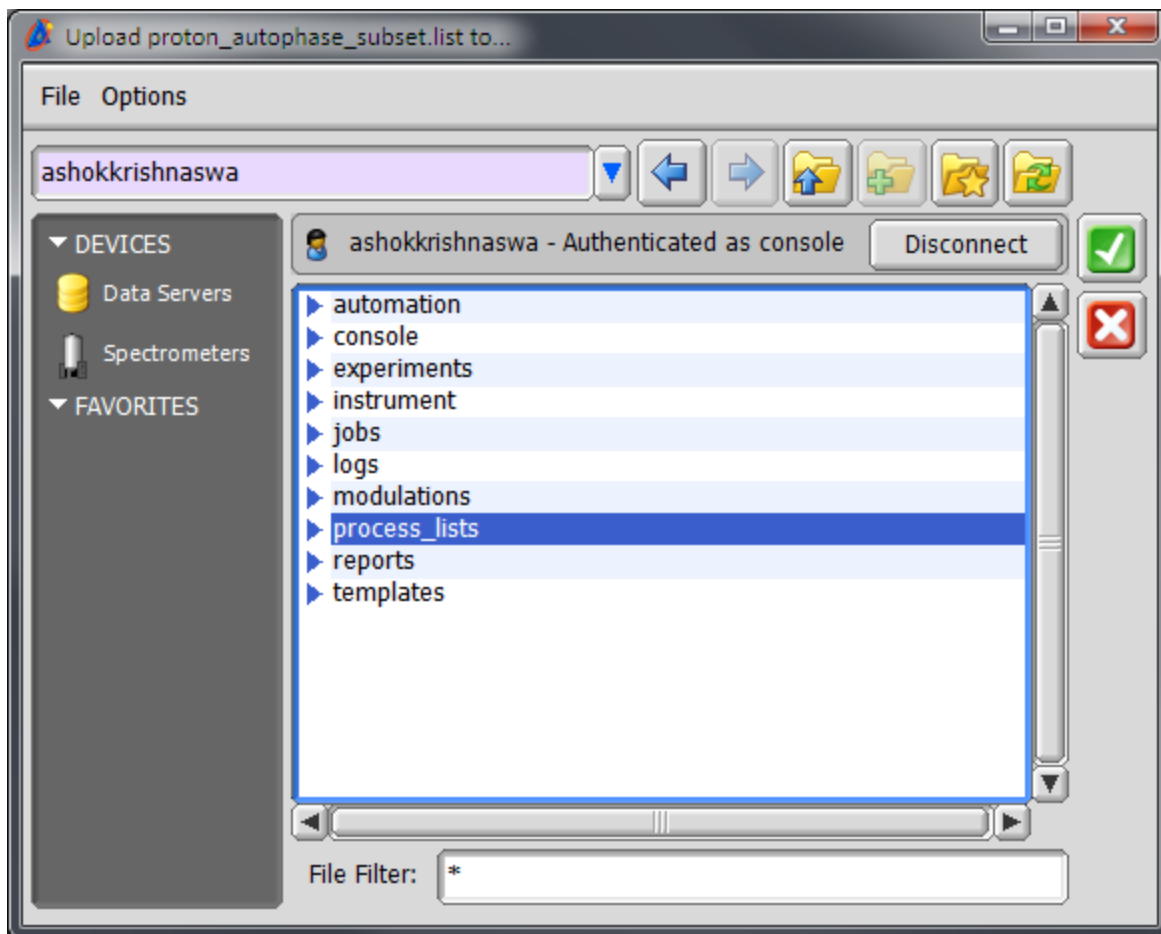


Double click on the "Process lists" directory.

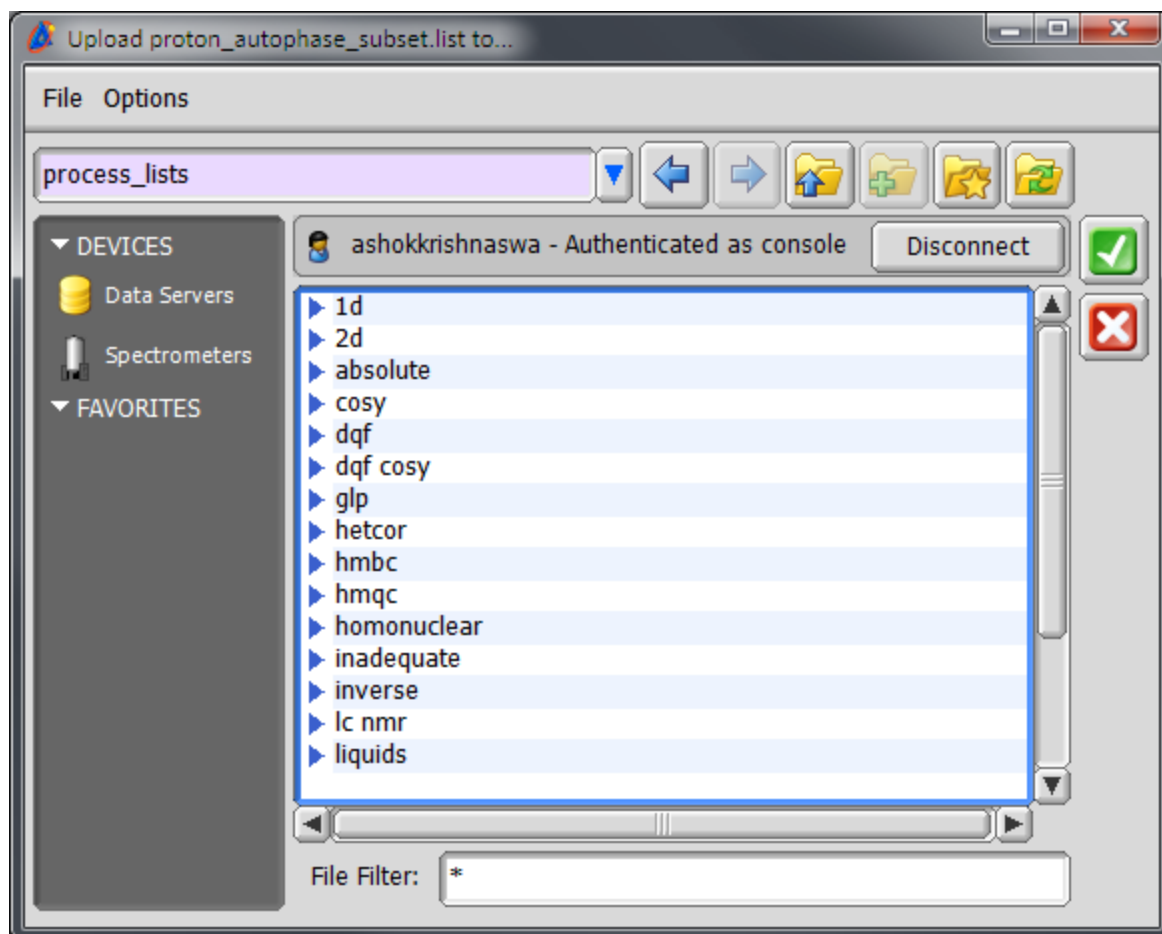

Hit the  button to upload the file.

The window open to the spectrometer will close.

Make sure that the message below (red outline) is displayed to confirm that the file uploaded successfully.

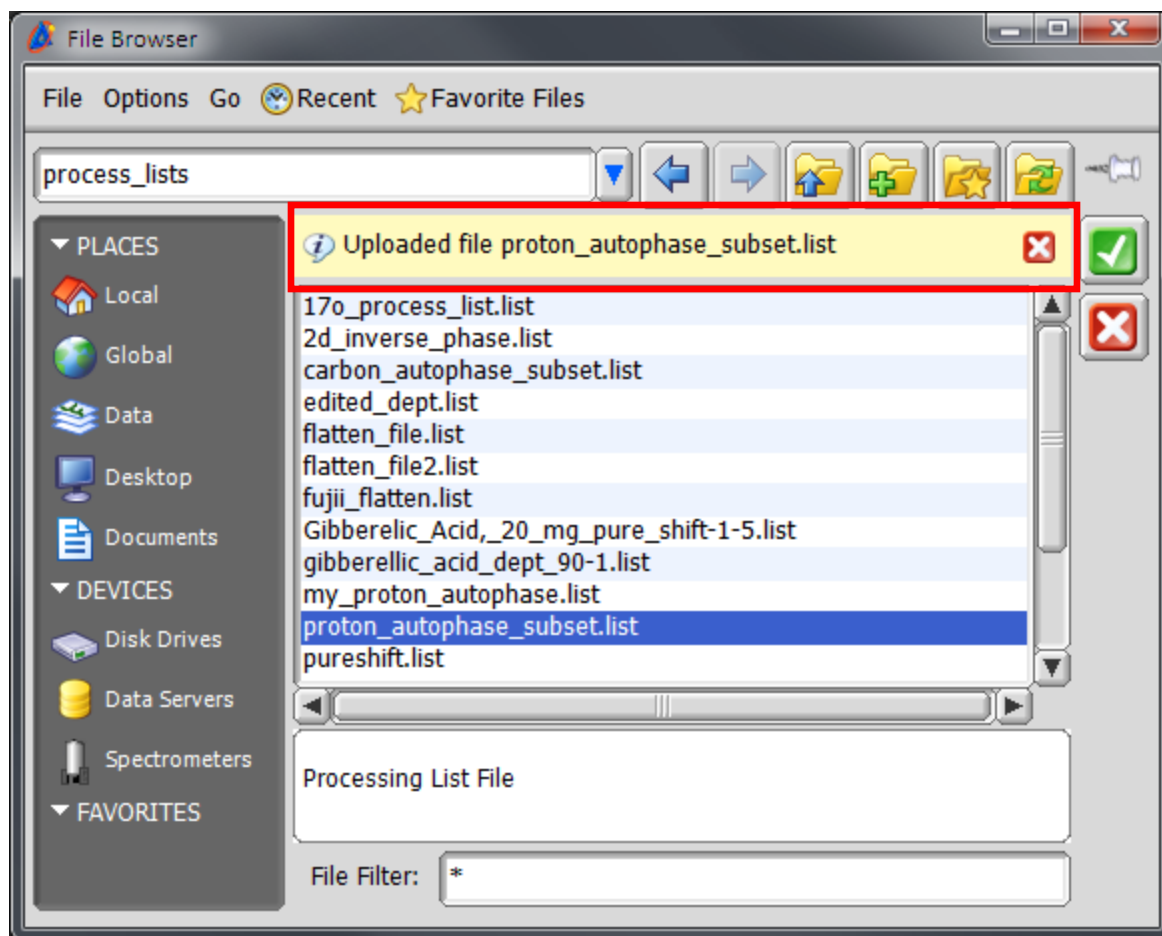

Repeat the process for “carbon_autophase_subset.list”.