



Pulling Log Files from a Cryomech CPA Helium Compressor Package

From the main menu screen, select the "Tools" menu

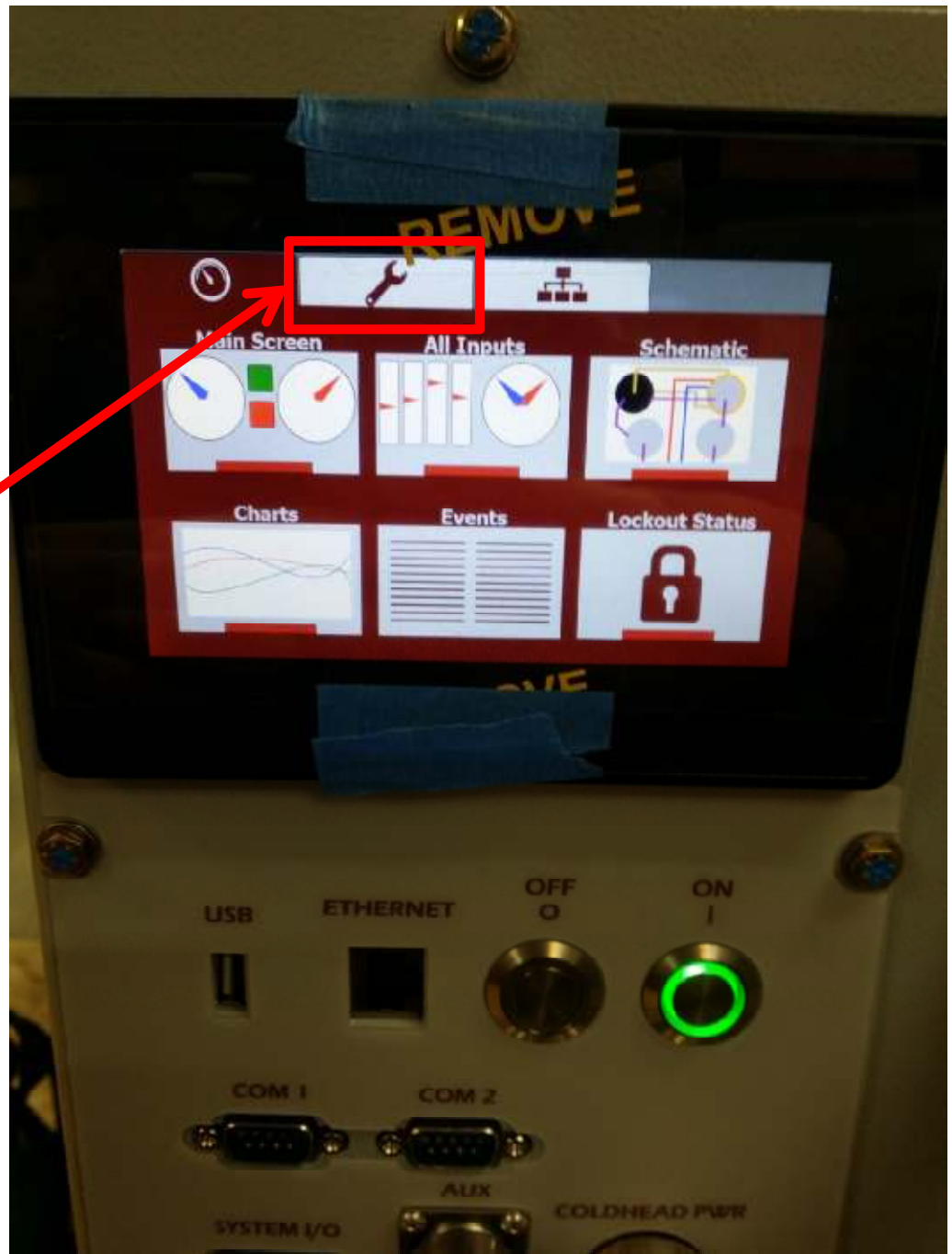

From the Tools menu, select the "Diagnostics" menu

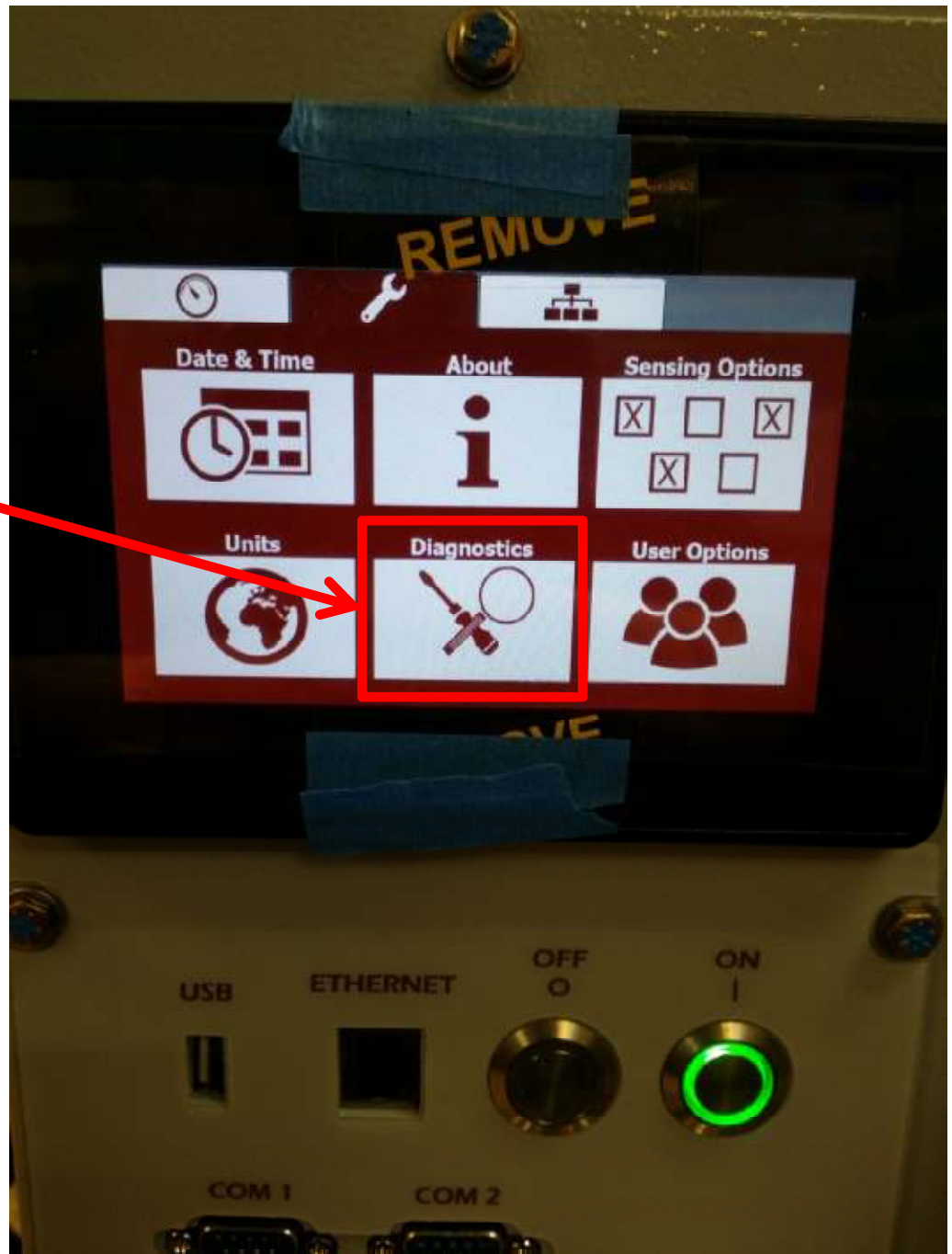

When in the diagnostics menu:

1. Insert Memory Stick into the USB port.
 - A USB with an LED indicator will be helpful for this.
2. Wait until the GREEN "USB device detected" button appears on the screen.
3. Once the button appears, select it and press OK to the next pop-up window.
4. Observe the LED light on the memory stick. Once it stops flashing, all of the files have been written to the device.
 - (4) .csv files will have been written to the device. These can then be emailed to Cryomech.
 - The device can then be removed from the CPA

