



Upload automation file to spectrometer


Disconnect from the Spectrometer before the following operation is performed.

Navigate to the location of the automation file directory. The extension for automation files is “JAF”. Select the automation file to be uploaded. In this example “Standard.jaf” is being uploaded.

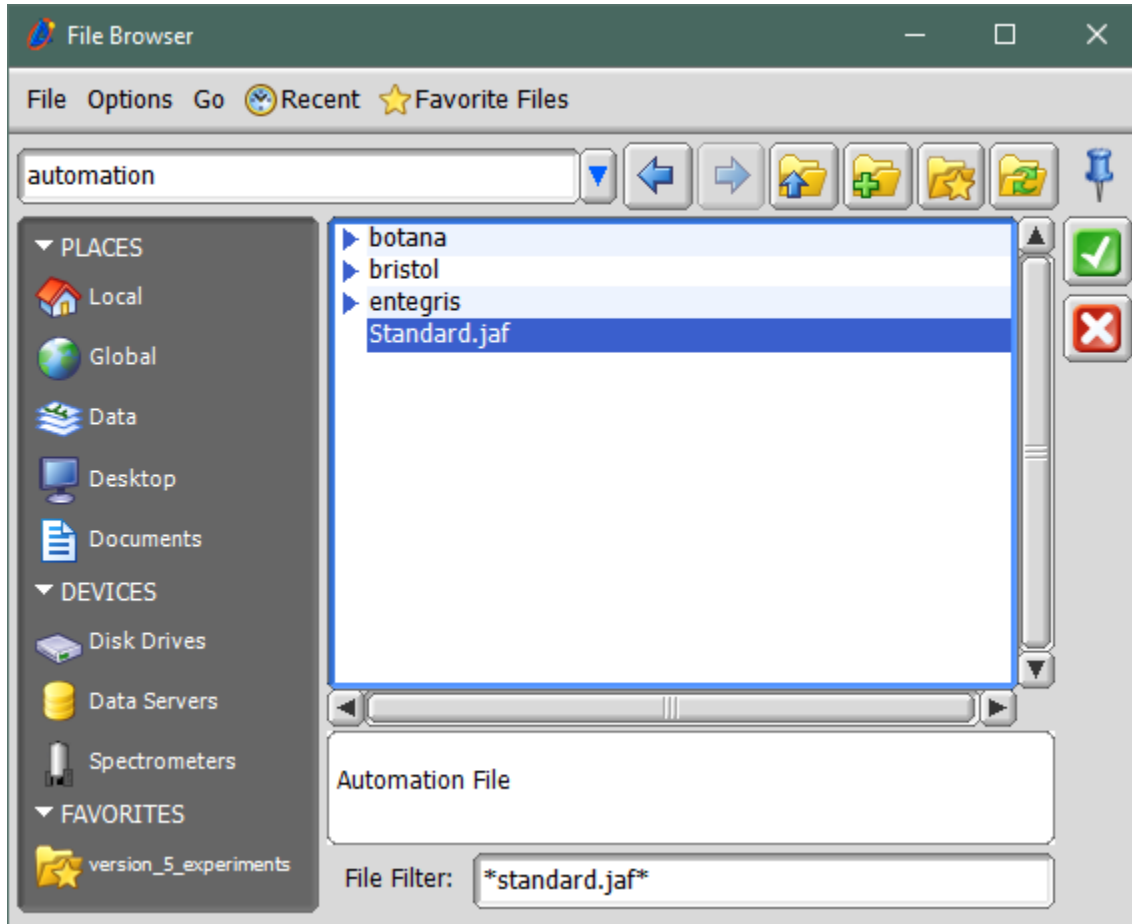


From the “File” menu select “Upload File”

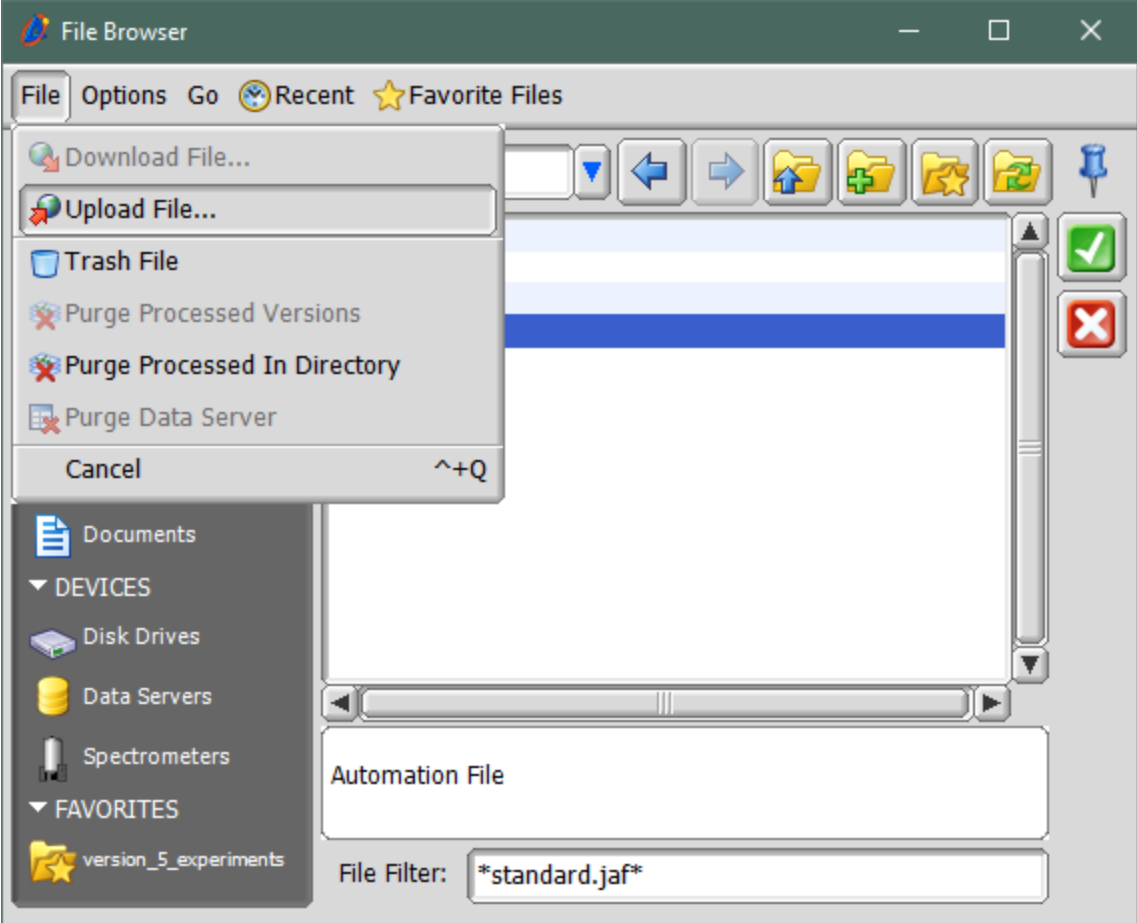


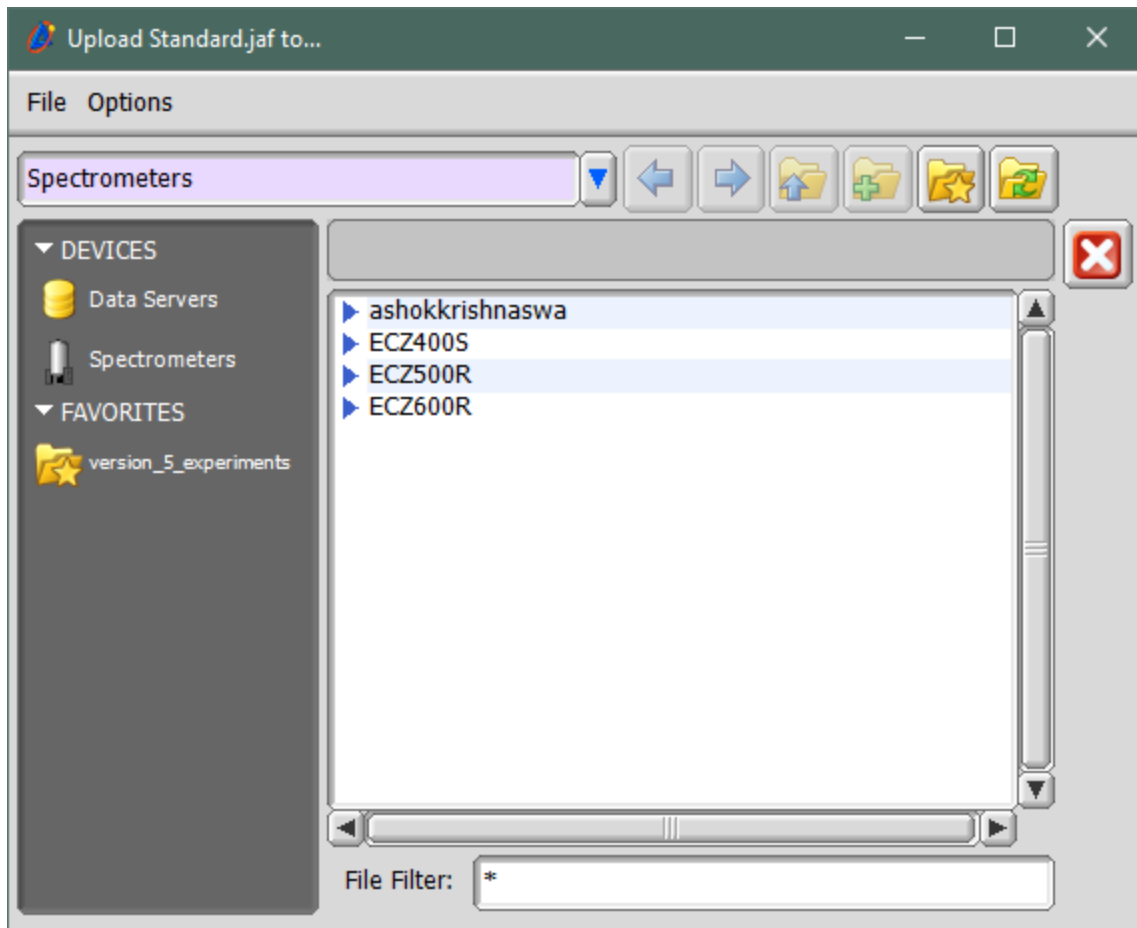

A new window displaying available spectrometers will appear. Select the spectrometer to which the file needs to be uploaded to.

Click on "Connect As..."

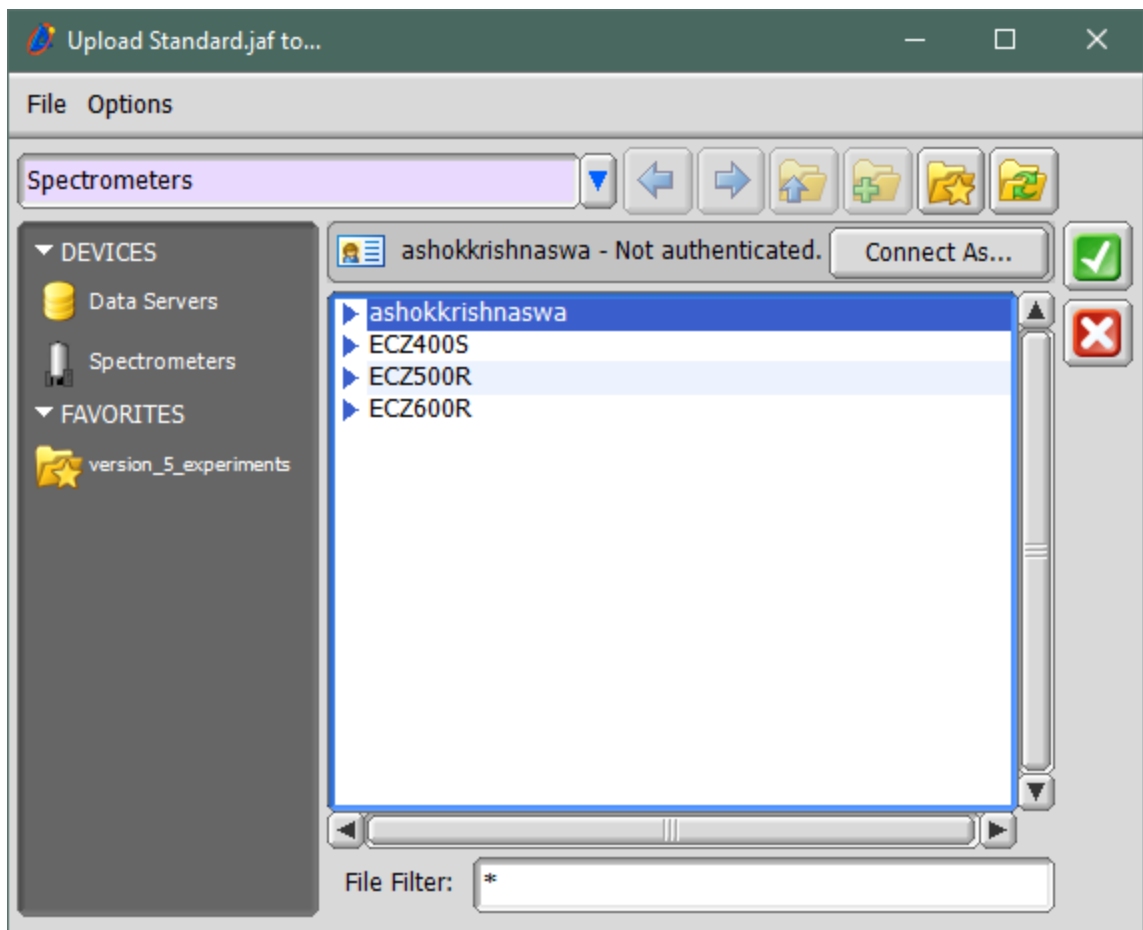

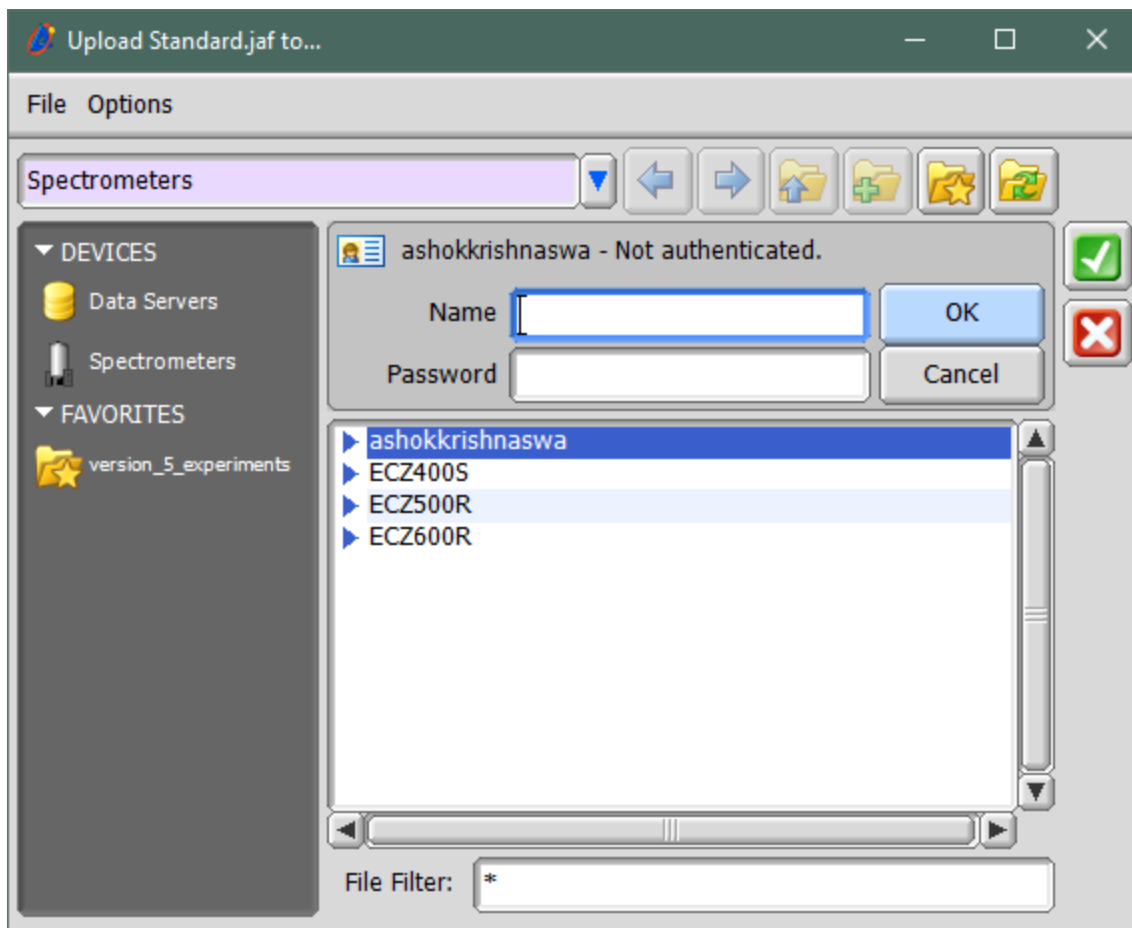


Enter


Name: console

Password: console (**Note: If the password has been changed, enter the current password for user “console”**).


Hit the “OK” button.

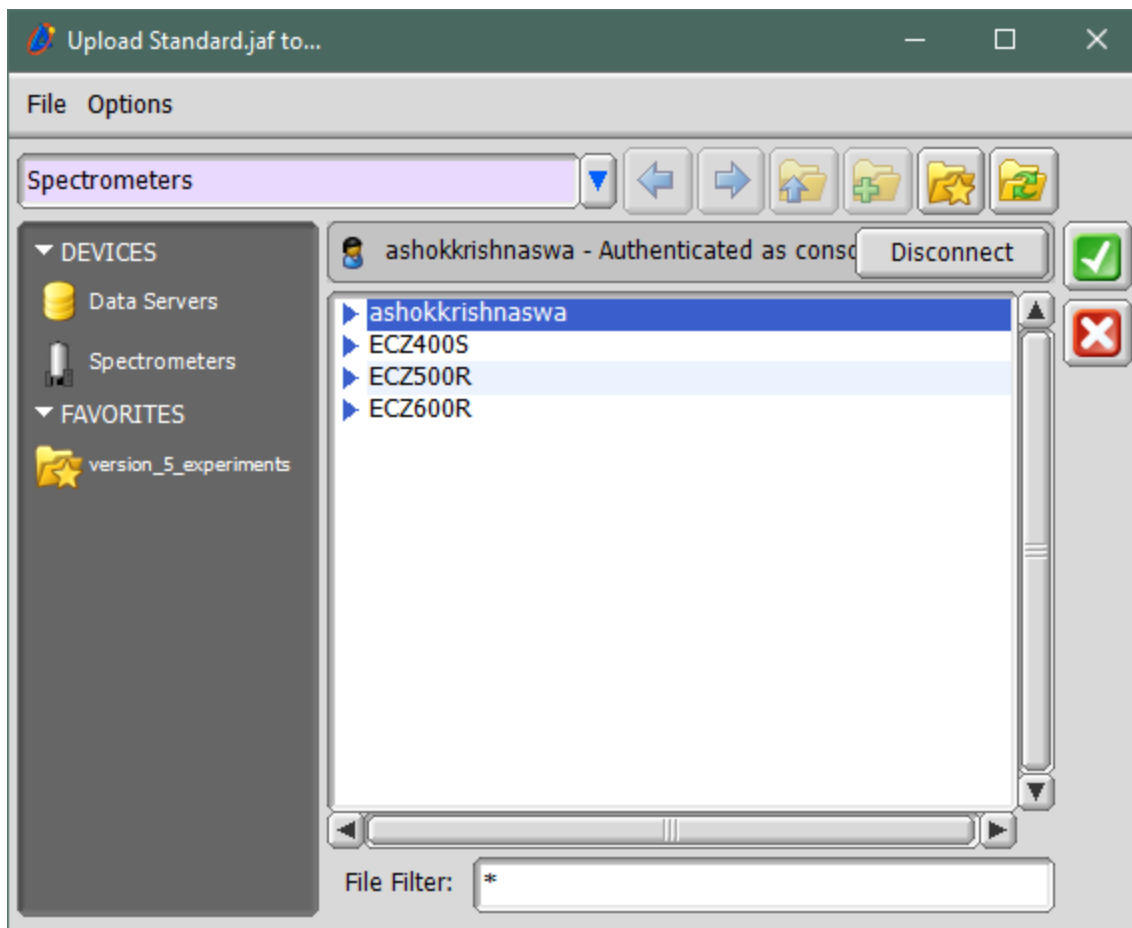


Double click on the selected spectrometer to display sub-folders.

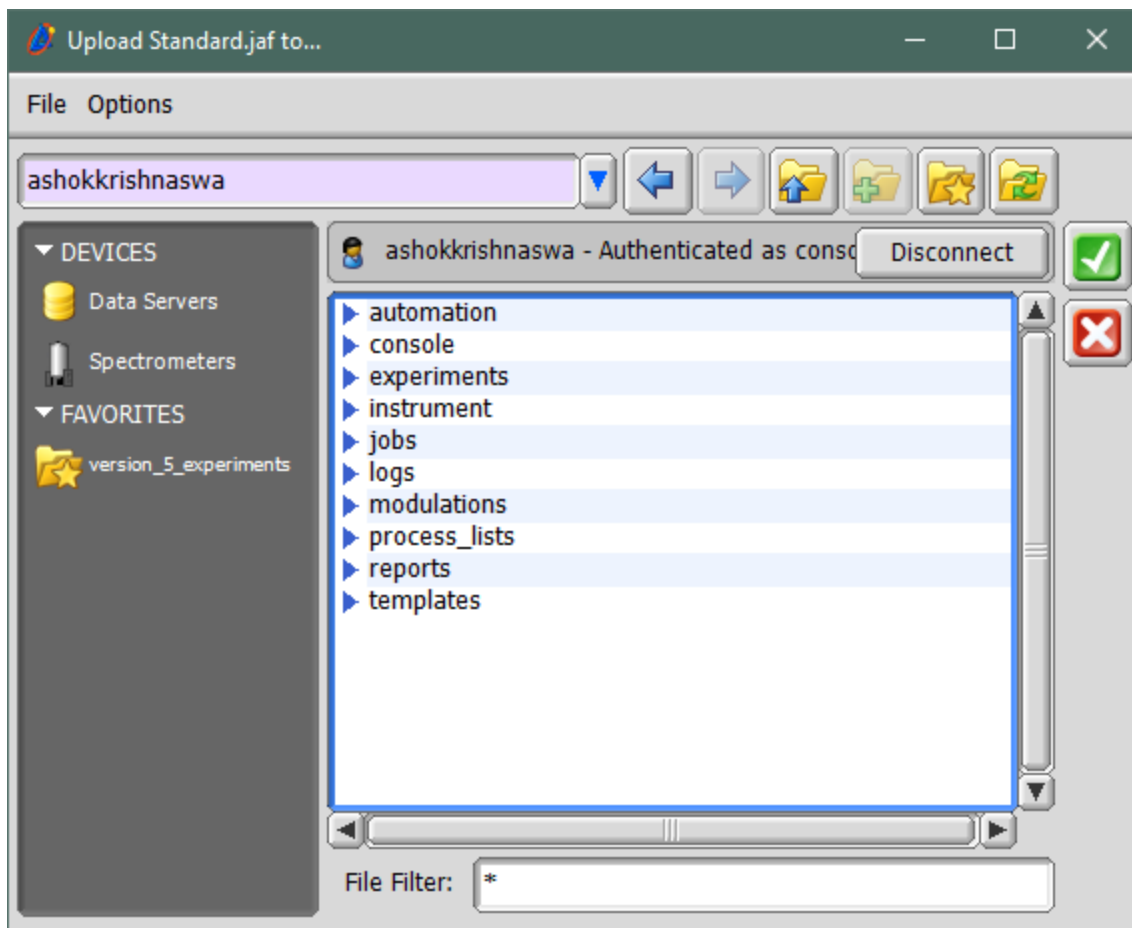

Double click on the “Automation” directory.

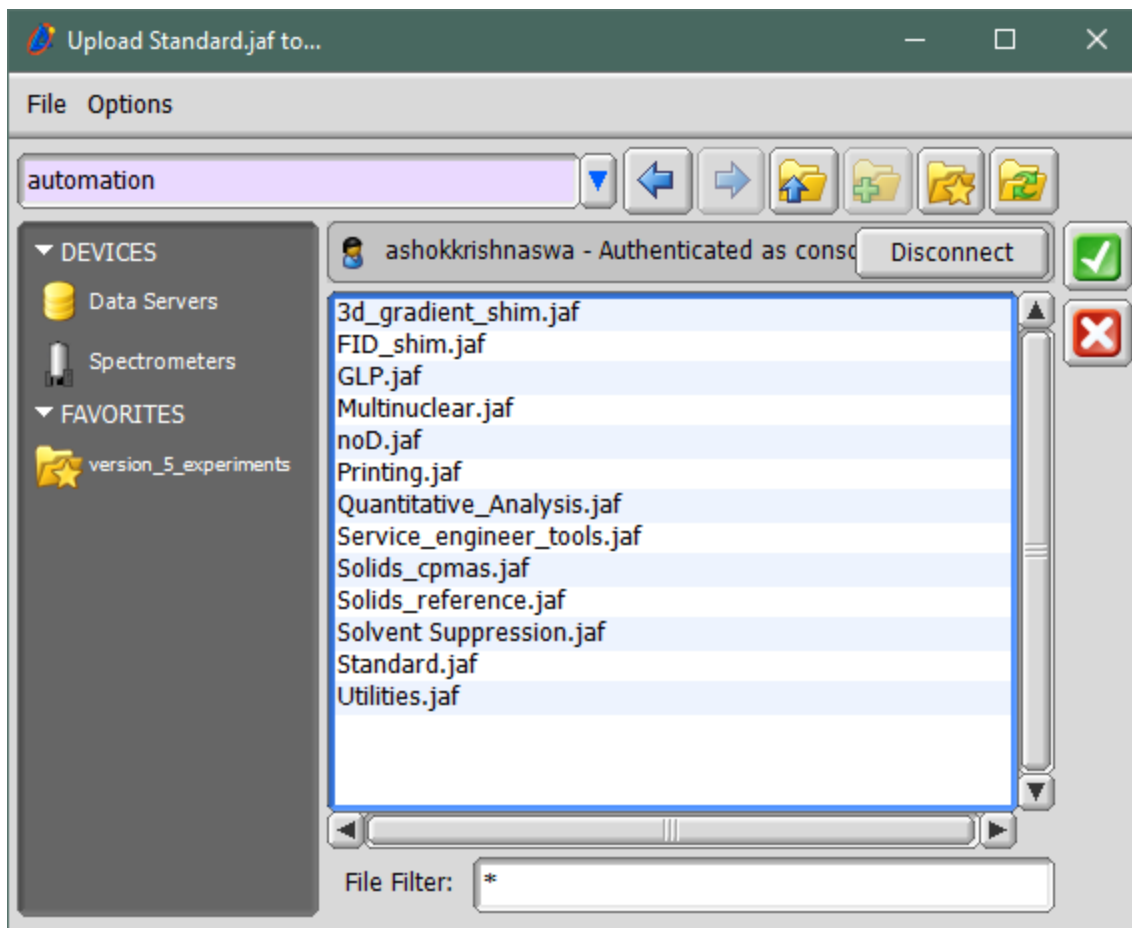

Hit the  button to upload the file.

The window open to the spectrometer will close.

Make sure that the message below (red outline) is displayed to confirm that the file uploaded successfully.

